



## ORBITAL AND LINEAR DIGITAL SHAKER

### USER MANUAL

CONTENTS	PAGE NUMBER
INTRODUCTION	2
WARRANTY	2
DELIVERY	2
SERVICE	2
1. SAFETY INSTRUCTIONS	2
2. PROPER USE	3
3. INSPECTIONS	3
4. CONTROL AND SET	4
5. OPEARTING MODES	5
6. SUPPORTED LOAD	7
7. FAULTS	8
8. INTERFACE	8
9. MAINTENANCE AND CLEANING	8
10. ASSOCIATED STANDARDS AND REGULATIONS	9
11. SPECIFICATIONS	10

## INTRODUCTION

Welcome to the Orbital and Linear Digital Shaker User Manual. Users should read this manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

PRODUCT CODE	DESCRIPTION
400.102.150	Linear Digital shaker (7.5Kg)
400.101.150	Orbital Digital shaker (7.5Kg)
400.102.100	Linear Digital shaker (2.5Kg)
400.101.100	Orbital Digital shaker (2.5Kg)

## WARRANTY

This instrument is warranted to be free from defects in materials and workmanship under normal use and service, for a period of 12 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of your local dealer.

## DELIVERY



Supplied with 1 x main unit, Power Cable and a user manual.

## SERVICE

You should contact your supplier for technical support.

- Please provide the representative with the following information:
- Product code
- Certification
- Description of problem (i.e., hardware or software)
- Methods and procedures adopted to resolve the problems
- Your contact information

## 1. SAFETY INSTRUCTIONS

	<p>Warning!</p> <p>Read the operating instructions completely before use and follow the safety instructions.</p> <p>Ensure that only trained staff work with the instrument.</p>
	<p>Protective ground contact!</p> <p>Make sure the socket must be earthed (protective ground contact) before use.</p>

Wear personal protective clothing/equipment in accordance with the hazard category of the media to be processed.

-Splashing of liquids.

-Vibration force resulting in broken glass containers/tubes.

- Use the instrument following the safety instructions and your own safety rules to avoid any accident.
- Users should not touch moving parts of the instrument to avoid any accidents that may occur. Pay attention to protect hands or fingers from rolling when moving the instrument.
- Set up the instrument in a spacious area on a stable, clean, non-slip, dry and fireproof surface. Do not operate the instrument in explosive atmospheres, with hazardous substances or under water.
- Observe the vessel on the shaking table for medium splashing out when setting the motor speed. Reduce the motor speed if the instrument is not running smoothly.
- Accessories must be securely attached to the instrument so that it cannot come off by itself and vessels must be firmly placed on the shaking table to avoid any injury.
- Working with combustible media or flammable materials may result in additional hazards. Only process media that will not react dangerously to the extra energy produced through using the instrument.
- Safe operation is only guaranteed with the accessories described in the "Accessories" chapter and by following the operating instructions in full. Always disconnect the plug before fitting accessories. Check the instrument and accessories beforehand for damage each time you use them. Do not use damaged components.
- The instrument may only be opened by trained technicians and always disconnect the plug beforehand.
- The operating voltage required as stated on the instrument and voltage in the supply network must match.
- Do not cover the instrument. Protect the instrument and accessories from bumps and impacts.
- Keep away from high magnetic fields.

## 2. PROPER USE

The instrument is designed for mixing or shaking liquids in schools, laboratories or factories. This instrument is not suitable for use in residential areas or other areas that may cause danger to the user or instrument as mentioned in chapter 1.

## 3. INSPECTION

### Receiving Inspection

Unpack the instrument carefully and check for any damages which may have arisen during transport. If this happens, please contact manufacturer for technical support.



**Note:** If there is any apparent damage to the system, please do not plug it into the power line.

### Trial Run

Do trial operation as follows:

- Make sure the required operating voltage and power supply voltage match.
- Ensure the socket is securely earthed.
- Plug in the power cable, Power ON the instrument.
- LCD displays the safe rotary speed limit.
- LCD displays operating mode.
- Turn the speed knob to set the rated rotary speed

- Press the speed knob and make sure the shaking function is switched ON.
- Press the motor regulation knob again and the shaking function is switched OFF.

If these operations above are normal, the instrument is ready to operate following the “operation” chapter. If these operations are not normal, the instrument may be in security protection state or be damaged.

If the mounting surface is not even, you can improve the stability of the instrument with the adjustable feet. To do this, turn the appropriate foot downward with the wrench until the instrument is standing securely on the surface.


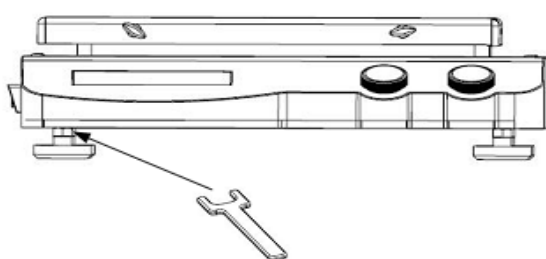


Figure 1

## 4. CONTROL AND SET

### Control

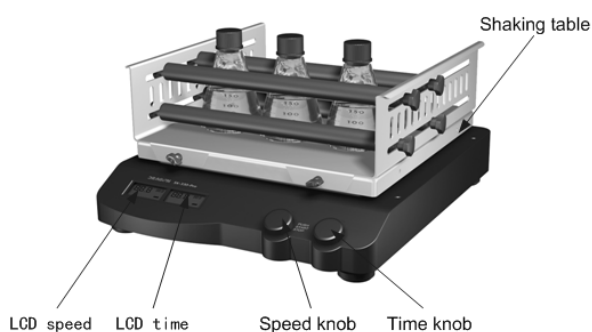

ITEMS	DESCRIPTIONS
<b>SPEED KNOB</b>	Set rated rotary speed, 'rotation' is switched ON or OFF via pressing the knob.
<b>TIME KNOB</b>	Set working time, "rotation" is switched ON or OFF via pressing the knob
<b>LCD</b>	Display the status of instrument and anything set
<b>MAINS SWITCH</b>	Switch ON or OFF.

Table 2

- Put the instrument on a stable and safe place and plug in the mains power.
- Turn ON the mains switch on the left panel.
- The instrument starts self-test.

- The instrument shows rated speed and time after initialisation.
- Turn the speed knob on the left side to set the rated speed.
- Press the speed knob and characters on the left LCD do not flash any more, and the shaking function is switched ON.
- Press the speed knob again and characters on the left LCD begin to flash, and the shaking function is switched OFF.
- Turn the time knob on the left side to set the rated time.
- Press the time knob and characters on the LCD do not flash any more, and the timing function is switched ON.
- Press the time knob again and characters on the right LCD begin to flash, and the timing function is switched OFF.

## Set time

The shaking time can be set by turning the time knob on the right. A distinction is made for the shaking time setting between timed mode and continuous operation. If continuous mode is selected, the shake can continue its shaking function for any amount of time with the previously set speed. If timed mode is selected, the shaker can run in the set time. After the instrument is powered OFF and restarted, the set time is erased to zero (blinking) and it switches into offline operating mode.

If a target time (max. 19h 59min) and speed are set, users can activate the instrument by pressing the time knob and the time to begin to read, if:

- Pressing the time knob, the speed and the time will stop. Pressing the time knob again, restart the shake function, the time will use the pre-set time.
- Pressing the speed knob, the speed and the time will stop; Pressing the time knob again, restart the shake function, the time not (There are flashing on the right of the LCD), pressing the time knob again restart the time function. The time will use the pre-set time.



**Note:** The current set time can be varied at any time.

## Set speed

The rated speed and upper speed limit can be set by turning the speed knob (As in Figure 6). The continuous mode can be switched ON with pressing the left knob without the time setting and switched OFF with pressing the knob again.

Adjust the motor speed knob slowly in order to keep the instrument running smoothly.



**Note:** The current set speed can be varied at any time.

## 5. OPERATING MODES

### Mode A

Operating mode A is the default mode set at the factory. After turning ON the instrument with the mains switch, the shaking and timing functions are turned OFF. The LCD displays the set time and speed values. They are adopted or varied when the corresponding functions are turned ON. After turning the instrument ON, the following process will be shown on the

display. The left LCD displays "SAF" and the right displays the upper speed limit (rpm) which can be set by pressing the speed knob and turning at the same time.



Figure 3

The LCD displays run modes "StA", "A" or "B" for about 2 seconds.



Figure 4

The set or stored speed value is shown in the left display, and the set or stored time value is shown in the right display. Turn the speed knob and time knob to set the speed and time.




Figure 5

After pressing the speed knob or time knob, the instrument begins to run at the set speed. The actual speed and time remaining are shown on the LCD. If the remaining has reached zero, the shaker stops its motion. Pressing the speed knob alone, the instrument will work in continuous mode (The time function is OFF)



Figure 6



**Note:** The set values can be varied during shaking. Shaking can be stopped by pressing the left or right knob.

## Mode B

When the instrument is turned ON and values for speed and time are erased to zero, users can set. The upper speed limit that has been set is taken over from operating mode A and cannot be modified. After the power has been turned OFF, the instrument will no longer automatically start up in operating Mode B by itself.

After turning ON the instrument and selected run mode, the following process will be shown on the display.

- The left LCD displays "SAF" and the right displays the upper speed limit (rpm) which cannot be changed.



Figure 7

- The LCD displays run modes "StA" for 2 seconds.



Figure 8

- The set speed value is shown in the left display, and the set time value is shown in the right display. Then set the motor rotary speed and time.



Figure 9



Figure 10

- After pressing the speed knob or time knob, the instrument begins to run at the set speed. The actual speed and time remaining are shown in the LCD. If the remaining has reached zero, the shaker stops its motion.



Figure 11



**Note:** The set values can be varied during shaking. Shaking can be stopped by pressing the left or right knob.

## Switching the mode

Switch the operating mode as follows:

- Turn the OFF button on the instrument with the mains switch.
- Hold down both knobs and turn ON the instrument with the mains switch. After 5 seconds, you can let go of the knobs.
- Switch between operating Modes A and B in order.

## 6. SUPPORTED LOAD

For safety reasons, the shaker must only be operated within the supported weight.

Make sure the holding surface of the shaker is always clean and level. Ensure the individual shaking vessels are to be placed evenly and fastened securely in the middle of the shaking table regardless of whether one or multiple shaking vessels are placed.



## 7. FAULTS

- When switched ON, the instrument doesn't work
  - Check whether the power cable is fitted well
  - Check whether the fuse is broken
- The speed cannot reach set value
  - Check whether it is overloaded
- The motor does not start via pressing the speed knob and time knob
  - Check whether the time is set to zero

## 8. INTERFACE

The Orbital and Linear Digital Shaker, special adapter and cable are used to link to external device and the standard 9-pin interface is to connect with the PC.

- The cable between the shaker and the computer uses EIA Standard RS232C communication line, corresponding to the DIN 66020 interface.
- Transmission method: Asynchronous signal transmission.
- Mode of transmission: Fully Duplex.
- 1 start bit; 8 character bits; 1 stop bit.
- Transmission speed: 9600 bit/s



**Note:** Do not plug the power cable in and out while the instrument is in use.

## 9. MAINTENANCE AND CLEANING

Use and maintain the instrument in correct way to keep it in a good operating state. This will lengthen its life time and avoid the damage to the instrument. Only use cleaning agents which have been approved by manufacturer to clean the instrument.

Only use cleanser as advised below:

<b>DYES</b>	Isopropyl alcohol
<b>CONSTRUCTION MATERIALS</b>	Water containing tenside / isopropyl alcohol
<b>COSMETICS</b>	Water containing tenside / isopropyl alcohol
<b>FOODSTUFFS</b>	Water containing tenside
<b>FUELS</b>	Water containing tenside

For materials which are not listed, please request information from manufacturer or your supplier. Before using anything other than the recommended method for cleaning or decontamination, the user must ascertain that this method does not destroy the instrument. Wear proper protective equipment during cleaning of the instrument.



**Note:**

- Electrical instrument may not be placed in the cleansing agent when purpose of cleaning.
- The instrument must be cleaned and put it into the initial packaging carton before sending for service or repair, avoiding the contamination of hazardous material.
- Switch OFF the instrument and put it in a dry, clean, stable place at room temperature when in long-term disuse.

## 10. ASSOCIATED STANDARDS AND REGULATIONS

**Construction in accordance with the following safety standards**

EN 61010-1

UL 3101-1

CAN/CSA C22.2(1010-1)

EN 61010-2-10

**Construction in accordance with the following EMC standards**

EN 61326-1

## 11. SPECIFICATIONS

ITEMS	SPECIFICATIONS
VOLTAGE [VAC]	100 ~240
FREQUENCY [HZ]	50/60
POWER [W]	30
SHAKING MOVEMENT	Orbital/ linear
ORBITAL DIAMETER [MM]	10 (7.5 kg) & 4.5 (2.5 kg)
MAX. SHAKING WEIGHT WITH ATTACHMENT	10kg (7.5 kg) & 4.5kg (2.5 kg)
MOTOR TYPE	External rotor brushless motor
MOTOR RATING INPUT [W]	28
MOTOR RATING OUTPUT [W]	15
PERMISSIBLE ON TIME [%]	100
SPEED RANGE [RPM]	100-350 (7.5 kg, Linear) & 100-350 (2.5 kg, Linear)
SPEED DISPLAY	100-500 (7.5 kg, Orbital) & 100-800 (2.5 kg, Orbital)
TIMER	LCD
TIMER DISPLAY	Yes
TIME SETTING RANGE [MIN]	LCD
RUN TYPE	1-1199
DIMENSIONS [D×W×H MM]	Time / Continuous operation
WEIGHT [KG]	420×370×100 (7.5 kg) 340×300×100 (2.5 kg)
PERMISSIBLE AMBIENT TEMPERATURE [°C]	13.5 (7.5 kg) & 8.1 (2.5 kg)
PERMISSIBLE RELATIVE HUMIDITY	5-40
PROTECTION CLASS ACC. TO DIN EN60529	IP21
RS232 INTERFACE	Yes

Table 3



Printed technical specifications are subject to change without notification  
Labco® is a registered trademark.

See the full product range online.

1800 052 226

[sales@labcoscientific.com.au](mailto:sales@labcoscientific.com.au)

[labcoscientific.com.au](http://labcoscientific.com.au)