

Allerflow | Gluten

[Learn more!](#) [Watch the Video](#)

Hit the play button!



AllerFlow Gluten is a rapid and convenient test for detection of gluten residue on surfaces as part of an allergen control program. Combining Hygiena's convenient sample collection swab design with classic lateral flow technology, AllerFlow Gluten makes gluten testing easier than ever before. Unlike other kits that contain several components, AllerFlow Gluten only consists of two parts – a sample collection device and a lateral flow cassette. The convenient sample collection swab device contains pre-measured extraction buffer for consistent results with minimal handling. After collection, the sample is poured directly into the lateral flow cassette fill well and results appear within 10 minutes.

Benefits

- Convenient and quick sample collection requires minimal training
- Pre-moistened swab enables reliable collection to ensure accurate results
- Simple two part kit includes all required materials - no pipettes, bottles, scales, or heating blocks required
- Pre-measured, self contained extraction buffer provides consistent results
- Easy-to-interpret lateral flow test results in 10 minutes
- Specific to the toxic fraction of gluten from wheat, barley, rye, and oat
- No cross-reactivity with soy, rice, or corn
- Up to 15 month shelf life at room or refrigerated temperatures (2-25°C) *
(* From date of manufacturing).

Sensitivity

- Sensitive to 5 µg/swab (5 ppm) gluten within 10 minutes

Positive Result



Negative Result



Part Number	Description	Quantity
HG-ALF-GL	AllerFlow Gluten: Gluten Surface Residue Test Kit Shop	25